Tamson Instruments Specification s

Specifications TV12LT

Tamson visibility bath 15 litres, ASTM D445 - IP 71 - ISO 3104 - IP EM PJ - IEC 61868



+	Small footprint
\oplus	Detachable front window
\oplus	Internal LED light
\oplus	Ultra-high stability
\oplus	RS232 communication
+	Bath drain
+	Built-in cooling -50°C / -58°F
+	4 places, small bath volume
+	Fast cool down to -20°C

Tamson visibility baths are specially designed for tests that require ultra-precise temperature control, or processes that need to be followed visually, e.g. viscometry, thermometer and sensor calibration, density and reaction rate measurement, etc. The window is heated preventing built up of condensate. The TV12LT is specially designed for kinematic viscosity determination of aviation fuels. Please see table $\mathring{3}$: "suggested set-up for determining the kinematic viscosity of aviation fuels". The TV12LT is able to cool down from ambient to -20°C (usual test temperature for aviation fuels) within 60 minutes.

Construction

The stainless steel construction ensures an exceptional stable bath temperature which is further improved by an ingenious stirring mechanism with baffle plates. All wetted parts are made of stainless steel and brass, providing resistance against all usual bath fluids. The cover of the bath has four round ø51 mm openings with lids, for suspending glass capillary viscometers in holders.

Item	Unit	TV12LT		
		-42+20°C -43.6+68°F	-42+80°C -43.6+176°F	-50*+20°C -58+68°F
P/N 230V/50Hz		00T0410	00T0425	00T0470
P/N 115V/60Hz		00T0415	00T0430	00T0475
P/N 230V/60Hz		00T0420	00T0435	00T0480
Reading	°C or °F	Menu selectable		
Interface		RS232		
Setting	[°C]	0.01		
Stability **	[°C]	stdev 0.004, min/max ± 0.014		
Uniformity **	[°C]	stdev 0.004, min/max ± 0.013		
Heating	[W]	500 +700		
Heaters		2		
Bath volume	[L]	15		
Cover		1 cover with 4 x ø51 mm openings		
Window	[mm]	255 x 230		
Opening bath	[mm]	250 x 98		
Depth	[mm]	300		
Length	[mm]	670		
Width	[mm]	425		
Height	[mm]	720		
Power	[Watt]	Nominal 800, Maximum 2100		
Ambient condition	[°C]	18 23		
CE All models conform to CE regulations				
* ± 2°C production variation B/N 00T0470, B/N 00T0475 and B/N 00T049				

± 2°C production variation P/N 00T0470. P/N 00T0475 and P/N 00T0480 Measured @-20°C in methanol

The bath is fitted with adjustable feet for leveling. Different covers can be used. The use of thermo insulated windows and window heating ensures clean sight. A permanent LED light is located in the top plate to supply clear light and guarantees optimal visibility inside the bath. A bath overflow outlet protects against expanding bath fluid or when the bath filling is too high.

Agitation

A vane type stirrer with maintenance free bearings moves the bath fluid past a special heater ensuring optimal control temperature and excellent uniformity. A baffle plate directs the fluid resulting in excellent temperature control and homogeneity.

Fine adjustment and offset

After the bath has become stable, the set point may be more accurately adjusted in the range of -5.00° to + 5.00°, if necessary. The absolute temperature can be adjusted with 0.005°C accuracy.

Conventional bath: Temperature stable within ±0.01°C after 40 minu

0:40

Time

[hrs:min]

Specifications TV12LT

Tamson visibility bath 15 litres, ASTM D445 - IP 71 - ISO 3104 - IP EM PJ - IEC 61868

-8.8°C

Femperature

-9.8°C

0:00

Span

Varies for different part numbers. Span lies from -50° C/ -58° F up to $+80^{\circ}$ C/176°F.

The bath conforms to CE-regulation. Further the bath is equipped with a mechanical over temperature device which trips when in case of malfunction the bath exceeds the pre-set maximum temperature. This feature guarantees safe around the clock operation.

Accuracy

Recovery from temperature dip

TV12LT bath

Conventional rectangle bath

Inside glass viscometer capillary


Accuracy @ -20°C

Methanol standard deviation ± 0.004°C min / max ± 0.013°C



Homogeneity


Methanol standard deviation ± 0.004°C min / max ± 0.013°C

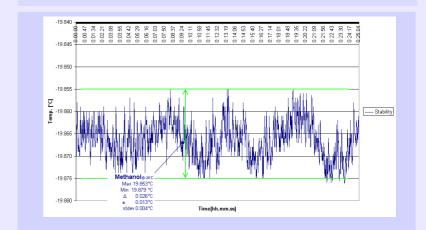

0:20

Cooldown


Reaches -20°C within 60 minutes Reaches -40°C within 120 minutes

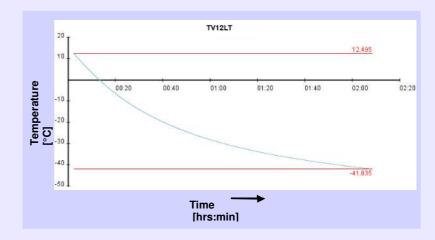


🔰 Contact: G-Labo Germany 吞 +49 6209 797100 🛕 info@g-labo.de 🏫 www.g-labo.de

Principle of operation

Superior homogeneity

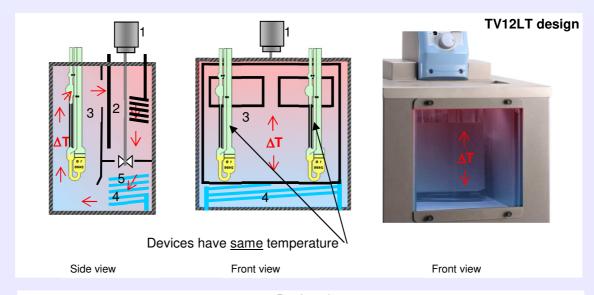

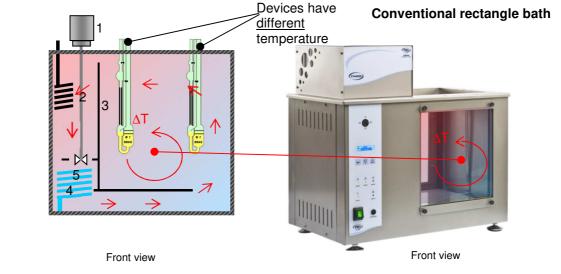


- 1: Motor
- 2: Heater
- 3 : Baffle plate
- 4 : Cooling
- 5 : Stirrer vane

Accessories

Table 1: TV12LT is standard included with:			
P/N	Picture	Description	
2272411		Cover with 4 openings: - 4 x ø51 mm opening - 2 x ø12.5mm opening for thermometer	
23T2411		4 * lid for ø 51 mm opening	

Table 2: Optional covers and levelling platform for TV12LT:		
P/N	Picture	Description
23T2413		Cover with 4 openings: - 4 x ø60 mm opening - 2 x ø12.5mm opening for thermometer
		4 * lid for ø60 mm opening
23T2412		Special cover for CFR (Cannon Fenske Routine) viscometers with 4 openings: - 4 x ø51 mm opening - 2 x ø12.5mm opening for thermometer
		4 * lid for ø 51 mm opening
13T6200		Levelling platform - without metal block (P/N 13T6210) - openings in cover can be custom designed (Please see page 6 for more information)
13T6210		Metal calibration block (Please see page 6 for more information)



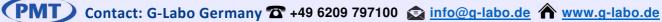

Tamson Instruments Specification sheet

Specifications TV12LT

Aviation Fuel Viscosity

Table 3: Suggested set-up to determine the kinematic viscosity of aviation fuel			
P/N	Picture	Suggested quantity	Description
00T0410	To the state of th		TV12LT (230V/50Hz)
00T0415		1	TV12LT (115V/60Hz)
00T0420			TV12LT (230V/60Hz)
10T6090	2.0 0 0 0 0 0 0 0 0 0 0 0 0 0 0 0 0 0 0	1	Timer, 8 positions
14T0303		1	Adapter to insert E20 thermometer in to the bath cover
19T4021		1	E20 thermometer Resolution 0.01 (two decimals) accuracy ± 0.02°C (Range -40°C to +140°C)
10T6001		4	Ubbelohde viscometer holder
25T0814	99	1	ISO 17025 Calibrated Ubbelohde viscometer tube size 0C
25T0816		1	ISO 17025 Calibrated Ubbelohde viscometer tube size 0B
25T0817		1	ISO 17025 Calibrated Ubbelohde viscometer tube size 1
25T0818		1	ISO 17025 Calibrated Ubbelohde viscometer tube size 1C
N2B		1	N2B viscosity reference standard
06T1724	171	1	Stoppers






Accessories

Table 4: Accessories and options			
Viscosity accessories		Please see specification sheet "Viscosity accessories", e.g. viscometers, viscometer holders, bath fluids, general purpose reference standards, etc	
02T0204		Spill tray Protects your lab against dripping and spilling during operation or when replacing bath fluid. The tray has a drainage valve 3/8" BSP connection.	
07T0085		Level detector/float	
13T3021		White contrast plate to get better visibility when measuring transparent liquids like jet fuel	





Accessories

Levelling platform (P/N 13T6200) and metal block (P/N 13T6210)

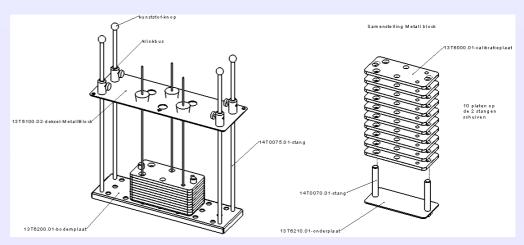

Dimensions top plate

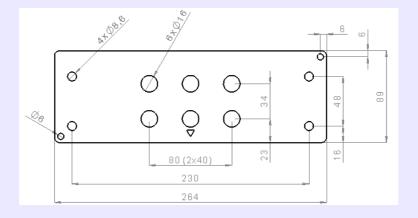

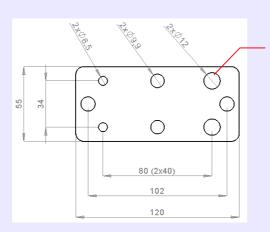

Number of holes, dimensions and position Other dimensions on request ®