Master Viscometers and Viscosity Oil Standards

ASTM D2162



- Ultra-stable ± 0.005°C
- Integrated cooling
- Stainless steel bath
- Heated window
- RS232 communication
- 0.001°C readout

General

This thermostatic bath is specially designed for tests that require ultra-precise temperature control, or processes that need to be followed visually. The TV16000 conforms to ASTM D2162: primary basic calibration of master viscometers and viscosity oil standards.

Construction

The stainless steel construction ensures an exceptional stable bath temperature which further is improved by an ingenious stirring mechanism with baffle plates. All wetted parts are made of stainless steel. This provides resistance against all usual bath fluids. The bath is fitted with four wheels, two of them having brakes. The cover of the bath is on request and has to be ordered separately. There's a wide range of possibilities such as round ø51 mm openings, ø61 mm openings, the use of a leveling platform or other square or round openings with other dimensions. A bath overflow outlet protects against overflow when bath fluid expands or raises due to submersion of parts.

Item	Unit	TV16000	
P/N 230V/50Hz		00T0490	
P/N 115V/60Hz		00T0492	
P/N 230V/60Hz		00T0495	
Power	[kW]	3	
Range	[°C]	10 60	
Reading		°C or °F	
Setting	[°C]	0.001	
Stability ±	[°C]	< 0.005	
Uniformity ±	[°C]	< 0.005	
Heating	[kW]	Control: 1 + Boost: 1.5	
Heating total	[kW]	2.5	
Bath volume	[L]	160	
Bath opening	[mm]	250 x 365	
Bath depth	[mm]	965	
Length	[mm]	575	
Width	[mm]	575	
Height	[mm]	1700	
Weight	[kg]	135	
Materials	Used inside bath: stainless steel 304, brass		
CE	Conforms to CE regulation		

Agitation

A vane type stirrer with maintenance free bearings moves the bath fluid past a special heater ensuring optimal bath temperature control and ensuring excellent uniformity inside the TV16000.

Span

Span lies from 10° C up to + 60° C (50° F / 140° F).

Safety

The bath conforms to CE regulation. Further the bath is equipped with a mechanical over temperature device which trips when in case of malfunction the bath exceeds the preset maximum temperature. This feature guarantees safe around the clock operation.

Optional

The system overall accuracy is within $\pm 0.005^{\circ}$ C based on measurement with water as bath fluid. This value is a min/max recorded over one hour. Standard deviation is better than $\pm 0.002^{\circ}$ C under these conditions.



Contact: G-Labo Germany 🛣 + 49 6209 797100 🙊 info@g-labo.de 🏠 www.g-labo.de

Master Viscometers and Viscosity Oil Standards

ASTM D2162

Fine adjustment and offset

After the temperature control is stable, the offset can be adjusted in the range of -5.00° C to $+5.00^{\circ}$ C, if necessary. The temperature can be set in 0.001°C resolution. A special offset offers ± 0.005°C resulting in setting of the set point temperature in ± 0.005°C.

Ambient condition

For proper cooling performance ambient temperature needs to be within the range of 18 .. 23°C.

Heating / cooling rate


Heating rate 12°C/hour Cooling rate 2.2°C/hour


Accuracy of control

Stability min / max over one hour measured in water, please see graphs on the next page @10 °C min / max ± 0.005 °C Standard deviation ± 0.0019 °C @20 °C min / max ± 0.003 °C Standard deviation ± 0.0013 °C @40 °C min / max ± 0.003 °C Standard deviation ± 0.0013 °C

Homogeneity ± 0.003 °C



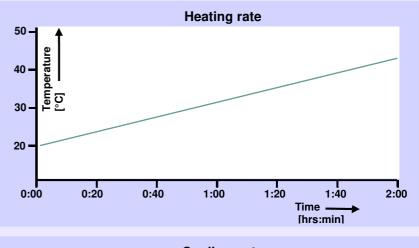


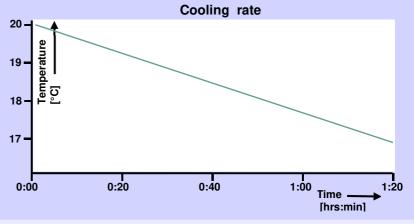





Master Viscometers and Viscosity Oil Standards

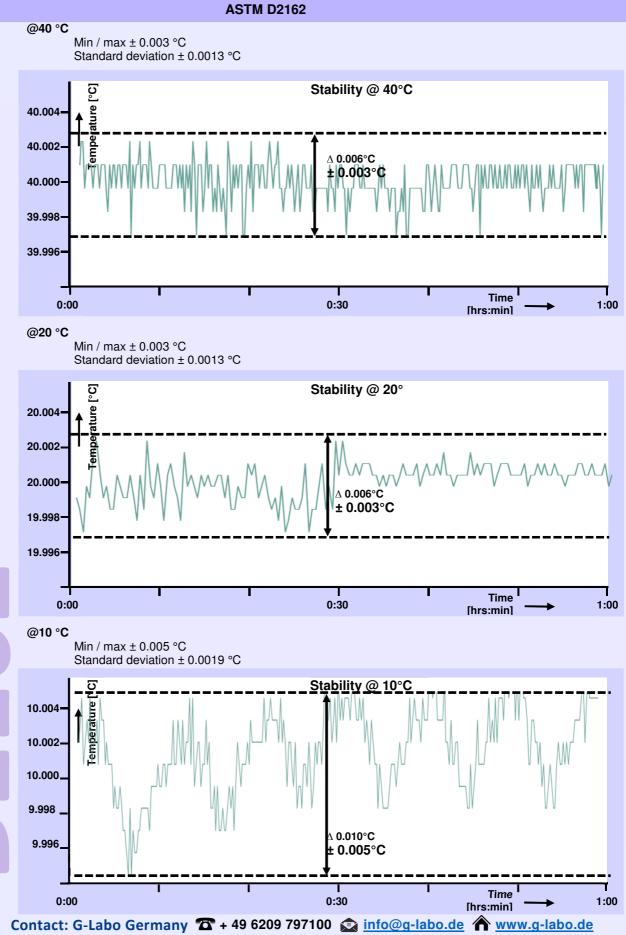

est 1878

Rev 1.1 - "17

Master Viscometers and Viscosity Oil Standards

ASTM D2162

	P/N	Picture	Description
	25T0581P	1	Master Ubbelohde viscometer ISO 17025 calibrated
-			under UKAS size 0 Master Ubbelohde viscometer ISO 17025 calibrated
	25T0583P	CP.	under UKAS size 0B
-	OFTOFOOD	20	Master Ubbelohde viscometer ISO 17025 calibrated
_	25T0582P	Q.	under UKAS size 0C
	25T0584P		Master Ubbelohde viscometer ISO 17025 calibrated
-			under UKAS size 1 Master Ubbelohde viscometer ISO 17025 calibrated
	25T0586P		under UKAS size 1B
-	OFTOFOED		Master Ubbelohde viscometer ISO 17025 calibrated
_	25T0585P		under UKAS size 1C
	25T0587P		Master Ubbelohde viscometer ISO 17025 calibrated
-			under UKAS size 2 Master Ubbelohde viscometer ISO 17025 calibrated
	25T0589P		under UKAS size 2B
-			Master Ubbelohde viscometer ISO 17025 calibrated
_	25T0588P		under UKAS size 2C
	25T0590P		Master Ubbelohde viscometer ISO 17025 calibrated
-			under UKAS size 3 Master Ubbelohde viscometer ISO 17025 calibrated
	25T0592P		under UKAS size 3B
-	25T0591P		Master Ubbelohde viscometer ISO 17025 calibrated
_	2010091P		under UKAS size 3C
	25T0593P		Master Ubbelohde viscometer ISO 17025 calibrated under UKAS size 4
-			Master Ubbelohde viscometer ISO 17025 calibrated
	25T0595P		under UKAS size 4B
	25T0594P		Master Ubbelohde viscometer ISO 17025 calibrated
-	20100041		under UKAS size 4C
_	25T0596P		Master Ubbelohde viscometer ISO 17025 calibrated under UKAS size 5
		1	
	10T6035		Viscometer holder Master Ubbelohde
	10T6030		Viscometer holder Master Cannon Fenske
		L	
	On request		Uncalibrated master viscometers either Cannon
			Fenske or Ubbelohde.





Master Viscometers and Viscosity Oil Standards

ASTM D2162

TV16000 accessories – options				
P/N	Picture	Description		
On request	~~~~	Top lid closed		
On request	000 000 000	Top lid ø51 mm openings		
On request	000 000	Top lid ø61 mm openings		

Alternative set-up TV7000DC + TLC15-5 with range from +15°C to +230°C and a temperature stability of \pm 0.007°C. For further information about the TV7000DC, please see specification sheet "TV7000DC".





PMT est. 1878

Contact: G-Labo Germany 🕿 + 49 6209 797100 🝙 info@g-labo.de 🏠 www.g-labo.de